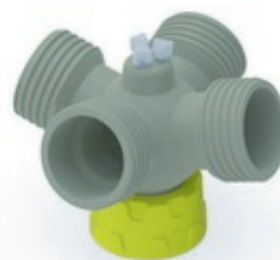


b.safe Distributor for Flushing Bottles Bohlender

[Product in eShop](#)

Home / [Brands](#) / [b.safe](#) / b.safe Distributor for Flushing Bottles Bohlender

Product Data

Screw cap made of PPS for GL45 thread / PP for GLS 80 thread, free movable insert made of PP with 4 lateral connectors (GL45) suitable for b.safe Caps GL 45 and 3 centric capillary connectors (UNF 1/4").

Applications

Flushing the HPLC installation without pollution or loss of solvents. The handling is very easy as the distributor is directly mounted on the flushing bottle. Remove the b.safe Caps from the eluent bottles and mount them on the connectors of the Distributor for Flushing Bottles. The passage at the distributor neck is sufficiently dimensioned so that the capillaries mounted on the cap can be inserted into the flushing bottle together with the filter. For the GL45 version of the Distributor for Flushing Bottles, we recommend to use smallest possible suction filters (max. O.D. 14 mm) to facilitate the insertion of tubing through the small bottle neck. If a connector of the distributor is not used, it can be closed with a screw cap GL45. Harmful solvent vapours cannot escape and the rinse solution is not polluted through ambient air by means of b.safe Caps with Air Valve. The Distributor for Flushing Bottles can also be used for back flushing the piston and/or the needle seat. Just connect the needed capillaries with b.safe Fittings on the centric capillary connectors as well as an Air Valve.

Item No.	for	Number of Blind Fittings	Number of flushing connectors	Number of capillary connectors	Price
94.M181.03	GL45	3	4x GL45	3	CHF 227.25
94.M181.10	GL80	3	4x GL45	3	CHF 262.10

More product images



* The prices are non-binding and are to be understood as selling prices in Swiss francs without value added tax (VAT), as well as all other fees, charges and taxes. The prices displayed in the eShop may differ from the PDF file due to regular updates.

** Please note that when ordering chemicals and detergents, transport and packaging costs for hazardous goods as well as legally prescribed fees are charged. These will be shown in detail on the order confirmation, which you will receive in addition to the confirmation of receipt.

*** Further information such as technical information and safety data sheets can be found online in our eShop.

**** The PDF file was created on www.huberlab.ch on 19.04.2026 at 17:07 o'clock.

www.huberlab.ch